GenCheck DNA Extraction kit [Type S/F]

For Soil and Fecal DNA



Product features

- Combined use of the chemical lysis method and the physical cell disruption method.
- Extracted from microorganisms with strong cell walls by Beads Beating.
- DNA can be extracted easily and quickly by a spin column.

Kit contents

1	Lysis Solution M	50 ml×1
2	Lysis Solution S	2 ml×1
3	Purification Solution	20 ml×1
4	Wash Solution 1	60 ml×1
(5)	Wash Solution 2	30 ml×1
6	TE(pH8.0)	5 ml×1
7	RNase A(100mg/mL)	100 μl×1
8	Beads Tube	50 pcs×1
9	Spin Column	50pcs×1





GenCheck DNA Extraction kit [Type S/F]

GenCheck® DNA Extraction Kit Type S/F is a kit for extracting DNA from soil and fecal samples. It combines the chemical lysis method with a surfactant and a physical cell disruption method by Beads Beating as the method to extract DNA. And, DNA is collected by centrifugation using a membrane based spin column used for collecting DNA. Therefore, DNA can be extracted easily and quickly.

Ex.1) Absorbance measurement result of DNA extracted with this kit

Sample	Place	N=3	Conc.(ng/µl)	λ260/λ280	λ260/λ230
	Tochigi, Japan	1	16.2	1.78	0.44
Andosol		2	15.2	1.67	0.44
		3	18.7	1.68	0.47
	Mt.Fuji, Japan	1	11.6	1.77	0.29
Volcanic ash soil		2	19.2	1.52	0.38
		3	16.2	1.6	0.34
	Nagasaki, Japan	1	35.4	1.8	1.11
Field soil A		2	40.9	1.84	1.4
		3	65.2	1.79	1.55
	Nagasaki, Japan	1	51.5	1.79	1.44
Field soil B		2	44.9	1.79	1.35
		3	41.4	1.78	1.09


Ex.2) Results of PCR using the extracted DNA of Example 1

- · ExTag polymerase/Takara
- ·27F/1492R
- Templete DNA: $lng/\mu l = 5\mu l = 20 \mu l = 20 cycle$

[Concentration measurement result by Qubit]

Sample	Place	N=3	Conc.(ng/µl)
	Tochigi, Japan	1	1.48
Andosol		2	1.95
		3	0.95
	Mt.Fuji, Japan	1	1.42
Volcanic ash soil		2	0.81
		3	2.04
	Nagasaki, Japan	1	2.99
Field soil A		2	2.79
		3	5.61
	Nagasaki, Japan	1	3.63
Field soil B		2	2.96
		3	4.03
		NC	Low

[Agarose electrophoresis(0.8%) *Load : 5 μ l]

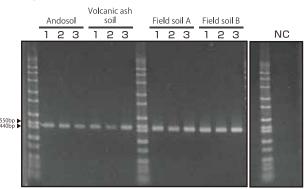

Product name	Catalog #	Size	Price
GenCheck® Extraction kit [Type S/F]	GC002-1N	50 prep	Contact

FASMAC CO.,LTD

5-1-3, Midorigaoka, Atsugi, Kanagawa, 243-0021 Japan

TEL: +81-46-295-8787 FAX: +81-46-294-3738

Web: http://fasmac.co.jp

